Experience perfection
through a combination
of quality, precision
and design!

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02. Liquid Handling · Dosiertechnik

No matter whether it is manual or automatic, aqueous solution or aggressive acid - we have the right device for your application!

Due to our long experience in the production of manual bottle top dispensers and automatic dosing systems, our products are easy to use, safe and precise. They are indispensable in everyday life in the laboratory.

We are proud to present our wide product range, which extends from bottle top dispensers and microliter pipettors to an electronic titrating device and an automatic dosing system. These cover most laboratory applications and a number of chemical and pharmaceutical manufacturing uses as well. With the widest product range on the laboratory market we offer real choice and expertise to our customers.

We can answer your dosing questions. Please contact us!

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bevorzugt empfohlen
- recommended

empfohlen
- not recommended

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* all media against which
polypropylene is resistant.

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For all other liquids please
contact us!

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FORTUNA® POLYFIX® Dispenser

FORTUNA® POLYFIX® - with glass piston

FORTUNA® POLYFIX® - with PTFE-coated piston

**FORTUNA® POLYFIX® - with amber glass cylinder
and glass piston**

**Due to the fact that PP is suitable for a lot
of chemicals, the POLYFIX® dispenser
can be used for most liquids.**

- visible dispensing procedure
- easy-to-use volume setting
system
- valve block made of
polypropylene with FIX adapter
- according to EN ISO 8655-5

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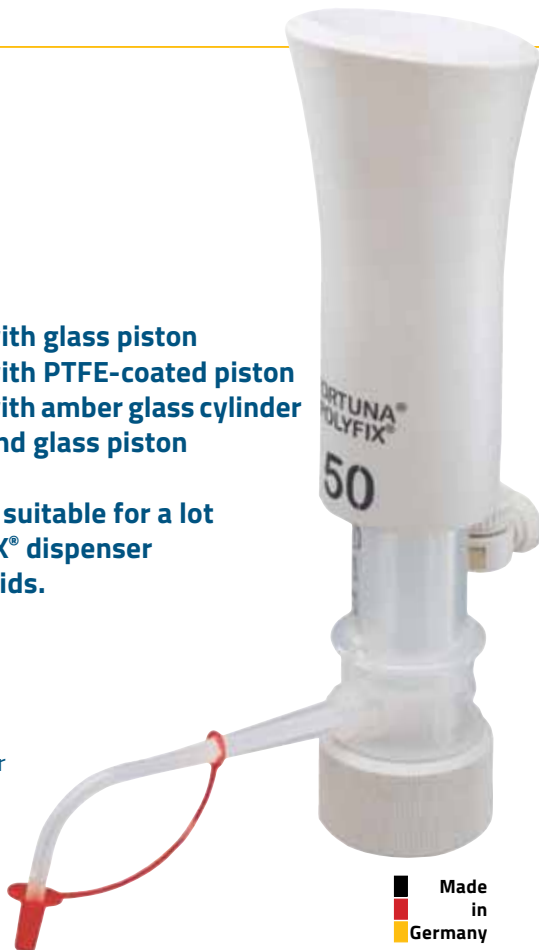
according to EN ISO 8655-5

101.070

- with glass piston
- suitable for all chemicals (incl. solvents), that do not attack polypropylene and don't crystallize

101.071

- with coated PTFE-piston
- suitable for acids and alkaline solutions that do not attack polypropylene

101.072

- with amber glass cylinder and glass piston
- suitable for all light sensitive liquids that do not attack polypropylene

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The bottle is not included in delivery (Matching bottles you find on p. 260 ff)

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comes with 1 PTFE
intake tube, 2 PP
adaptors (GL 40 + 45),
packing unit: 1

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Replacement Parts - for 101.070

101.077

Glass piston, complete with finger guard and volume setting system

101.075

Valve block unit, complete with dispensing cylinder and discharge tube

101.089

Stopper caps for discharge tubes, natural

101.093

Intake tubes, made of PTFE

- for 101.071

101.078

PTFE coated piston, complete with finger guard and volume setting system

101.075

Valve block unit, complete with dispensing cylinder and discharge tube

101.089

Stopper caps for discharge tubes, natural, pack 10

101.093

Intake tubes made of PTFE

- for 101.072

101.077

Glass piston, complete with finger guard and volume setting system

101.079

Valve block unit, complete with amber dispensing cylinder and discharge tube

101.089

Stopper caps for discharge tubes, natural

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101.093

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Replacement Parts for POLYFIX® Dispenser

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FORTUNA® OPTIFIX® Dispenser

Advantages at a glance

- **Best safety by use of high grade material**
- **Visible dosing procedure**
- **No contamination by metal parts, e.g. springs**
- **High accuracy, according to EN ISO 8655-5**
- **Premium quality**
- **BASIC and SOLVENT - wide volume range up to 500 ml**

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02. Liquid Handling · Dosiertechnik

Dispensing Piston

- proven technology
- made of glass, solid PTFE-encased, exception: SOLVENT
- virtually indestructible

Precision Dispensing Cylinder

- made of glass, dosing procedure visible (except type HF, made of PTFE)
- coated with plastic - protection against breakage
 - graduation clear readable
 - easy to dismantle and to clean

Valve Block

- 100 % tightness
- no metal springs - no contamination of the dosing liquid

Finger Protection Cap

- ergonomic type
- timeless design
- with colour-code-system

Volume Setting System

- fast and easy volume setting
- additional fine adjustment
- for right- and left hand use
- easy to read scale

FIX Adaptor

- to align the dispenser on a bottle
- made of PP for BASIC + SOLVENT
- made of PTFE for SAFETY +HF

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according to EN ISO 8655-5

101.080

**for dispensing light acids,
bases, aqueous solutions
and crystallizing liquids**

- with solid PTFE-encased dispensing piston
- dispensing cylinder made of glass
- visible dosing procedure
- no metal spring, no contamination of the dosing liquid
- volume up to 500 ml

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101.080
200 - 500 ml

**comes with 1 PTFE
intake tube, 2 PP
adaptors (GL 40 + 45),
packing unit: 1**

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101.081

for dispensing solvents

- with glass piston
- dispensing cylinder made of glass
- visible dosing procedure
- no metal spring, no contamination of the dosing liquid
- volume up to 500 ml

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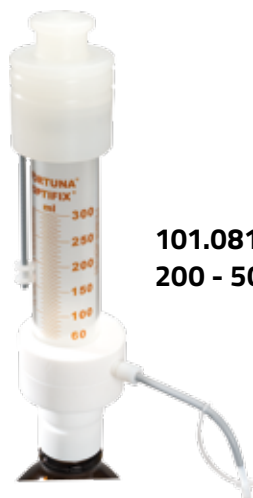
according to EN ISO 8655-5

101.106

**For all liquids that can
be handled with the
OPTIFIX dispensers
and for hydrofluoric acid**

- made of PTFE
- with PTFE-encased dispensing piston and PTFE-cylinder
- FIX adaptor made of solid PTFE
- no metal spring, no contamination of the dosing liquid
- volume up to 50 ml

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The bottle is not included in delivery (Matching bottles you find on p. 274)

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comes with 1 PTFE intake tube, 2 PTFE adaptors (GL 40 + 45), packing unit: 1

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according to EN ISO 8655-5

101.107

for dispensing strong acids and aggressive chemicals

- PTFE-encased dispensing piston with sealing lip
- dispensing cylinder made of glass
- visible dosing procedure
- FIX adaptor made of solid PTFE
- no metal spring, no contamination of the dosing liquid
- volume up to 100 ml

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comes with 1 PTFE
intake tube, 2 PTFE
adaptors (GL 40 + 45),
packing unit: 1

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according to EN ISO 8655-5

101.107 S

for dispensing strong fuming acids and very abrasive liquids

- with safety stopcock and additional safety valve in the ventilation tube
- PTFE-encased dispensing piston with sealing lip
- dispensing cylinder made of glass
- visible dosing procedure
- FIX adaptor made of solid PTFE
- no metal spring, no contamination of the dosing liquid
- volume up to 100 ml

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- for 101.080

101.000

Piston BASIC, solid PTFE encased, with finger guard and volume setting system

101.004

Valve block unit with dispensing cylinder

- for 101.081

101.001

Piston SOLVENT, made of glass, with finger guard and volume setting system

101.004

Valve block unit with dispensing cylinder

- for 101.107

101.003

Piston SAFETY, solid PTFE encased, with sealing lip, finger guard and volume setting system

101.004

Valve block unit with dispensing cylinder

- for 101.107-..S

101.003

Piston SAFETY S, solid PTFE encased, with sealing lip, finger guard and volume setting system

101.004-S

Valve block unit with dispensing cylinder, with safety valve

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Replacement Parts for OPTIFIX® Dispenser

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Volume setting system

101.000-KS

Volume setting system for BASIC, SOLVENT, HF, SAFETY, SAFETY S, 2-100 ml

101.000-KM

Volume setting system for BASIC, SOLVENT, 200 - 500 ml

Volumen-Einstellsystem

101.000-KS

Volumen-Einstellsystem für BASIC, SOLVENT, HF, SAFETY, SAFETY S, 2-100 ml

101.000-KM

Volumen-Einstellsystem für BASIC, SOLVENT, 200 - 500 ml

Intake tubes

101.093-36

Intake tubes for HF, 5 ml, made of PTFE

101.093-38/-48/-51/-58

Intake tubes for BASIC, SOLVENT, HF (10+30 ml), SAFETY, SAFETY S, made of PTFE

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10.1433

Tubing, made of PTFE, sold in varying lengths (state length with order)

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Discharge tube sets

101.089

Stopper caps for discharge tubes,
natural

101.094

Discharge tube set (tube with stabiliser,
fixing screw, cap)

101.095

Discharge tube set with Safety-Stopcock
(tube with stabiliser, 2 fixing screws,
PTFE-stopcock with dispensing tip),
completely interchangeable;
prevents dropping after the dispensing
procedure

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aus PTFE

101.105

Humidity absorber, made of PTFE, for hygroscopic liquids, single pack

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101.104

**for OPTIFIX® SAFETY,
SAFETY S, OPTIMAT®**

- flexible remote dispensing
- suitable for easy serial dispensing
- tubing and dispensing valve made of PTFE
- with new metal spring, chemical resistant
- prevents liquid from coming out and stops liquid flowing back

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**The WANDERER
for flexible
remote dispensing**

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FORTUNA® OPTIFIX® titrier

Benefits

- Easy to use
- easy to read LCD-display
- visible dispensing process
- left or right hand operation
- no metal spring
- high precision due to high tech micro electronic system
- can be used to dispense aggressive liquids

Application

- in chemical laboratories
- in environmental laboratories
- in food laboratories
- for routine daily work

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Technical Features

- with certification of conformity
- GS = safety approved
- cylinder cap. 50 ml, display max. 99.99 ml
- subdivision 0.01 ml, Accuracy = $\pm 0.6\%$, Tolerance = $\pm 0.2\%$
- according to EN ISO 8655-5

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LCD-display and buttons

- the size of the LCD-display figures is 10 mm
- the arrow direction indicators show clearly the FILL or TITration mode
- LCD-display and buttons are at 10° angle to ensure easy reading and operation
- the mode is switched automatically from FILL to TITration

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The dispensing unit

- the high grade registered glass cylinder is amber stained to allow use with light sensitive materials
- dispensing process is always visible and, if necessary, air bubbles can be easily removed
- the precision dispensing piston, made of glass, is coated with solid PTFE
- excellent resistance both to aggressive acids and alkaline solutions

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The valve block system

- made of pure PTFE
- can be directly mounted on bottles with GL 32
- has an additional ball valve in the ventilation tube

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according to EN ISO 8655-5

101.102

**for dispensing strong acids,
alkalis and aggressive liquids**

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101.102EEZ

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101.094-48T

**Discharge tube ,titrier', with
extra thin tip for titrating drop
by drop**

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FORTUNA® OPTIMAT®
dispensing pump

101.126

for dispensing strong fuming acids and very abrasive liquids, single pack (intake and discharge tube sets are not included)

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Pumps for OPTIMAT® programmable dispenser:

Packing unit: 1

Pumpen für OPTIMAT® Dosierstation:

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					101.126
Volume	Grad.	A* ≤ ±	CV* ≤ ±	max.V	Art.-No.
Volumen	Grad.	R* ≤ ±	V* ≤ ±	max.V	Art.-Nr.
0 - 1 ml	0.05 ml	± 0.6 %	± 0.2 %	0,55 ml/s	101.126-20
0 - 2 ml	0.1 ml	± 0.6 %	± 0.2 %	1,90 ml/s	101.126-27
0 - 10 ml	0.2 ml	± 0.6 %	± 0.2 %	3,35 ml/s	101.126-38
0 - 30 ml	0.5 ml	± 0.6 %	± 0.2 %	9,60 ml/s	101.126-44
0 - 50 ml	1.0 ml	± 0.6 %	± 0.2 %	15,20 ml/s	101.126-48
0-100 ml	2.0 ml	± 0.6 %	± 0.2 %	24,30 ml/s	101.126-51
0-200 ml ⁽¹⁾	5.0 ml	± 0.6 %	± 0.2 %	31,30 ml/s	101.126-53
0-300 ml ⁽¹⁾	5.0 ml	± 0.6 %	± 0.2 %	31,80 ml/s	101.126-56

* A=Accuracy, CV=Coefficient of variation *R=Richtigkeit, V= Variationskoeffizient

⁽¹⁾ additional to ISO 8655 · in Ergänzung zur ISO 8655

02. Liquid Handling · Dosiertechnik

Accessories for OPTIMAT® and OPTIMAT® dosing pumps

101.132
Intake tube
length 1 m

101.134
Discharge tube set
made of PTFE, length 0.75
or 1.50 m

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155.000-40
Serielles Verbindungskabel

**Tubes for
150.000-3 OPTIMAT®
programmable dispenser:**

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